

INSTRUCTIONS FOR USE

ZEO-Z009



CAUTION: Laboratory spills may contain broken glass. Avoid direct contact and use the supplied pan and scraper to reduce the risk of cuts.

SAFE SPILL RESPONSE PROCEDURE

Complete each step in sequence

1 CONTAIN THE SPILL

Apply **Absorbent Powder** around the outside perimeter of the spill first to contain the fluid. Sprinkle generously over the remainder of the spill and allow approximately **3-5 minutes** to congeal.

2 APPLY PERSONAL PROTECTIVE EQUIPMENT

Don the relevant **gloves, protective eyewear, mask and plastic apron**.

3 COLLECT BROKEN GLASS SAFELY

Laboratory spills may contain broken glass. Avoid direct contact and cuts. Use the supplied pan and scraper to collect broken glass, then place it into the plastic container (**10 L bucket**).

4 COLLECT THE ABSORBED MATERIAL

Using the disposable plastic pan and scraper, collect the absorbed substance and place it into a **yellow contaminated-waste bag**. Secure the bag tightly using one of the supplied ties.

5 REMOVE RESIDUAL CONTAMINANT

Wipe the affected area with absorbent cloths to remove any remaining contaminant.

6 CONTAIN ALL USED ITEMS

Place all used items in a yellow waste bag and seal it. The original **10 L bucket**, with the lid securely fitted, may also be used to contain spill waste, particularly where broken glass is present.

7 DISPOSE IN ACCORDANCE WITH REGULATIONS

Dispose of contaminated product in accordance with applicable **State/Territory regulations**.

8 CLEAN THE SPILL AREA

Wash the spill area with warm, soapy water.

9 COMPLETE HAND HYGIENE

Wash hands thoroughly after cleaning is complete.

10 REORDER THE REPLACEMENT KIT

Reorder a replacement kit after use. **Reorder code: ZEO-Z009.**



Completion check: Confirm contaminated materials are sealed for disposal, the spill area has been washed, hand hygiene is complete and a replacement kit has been ordered.

V 202605 - The images and/or information provided in this document are for general guidance only and may be subject to technical changes. We reserve the right to modify product specifications, design, or information where necessary. Images are for illustrative purposes only and may differ from the actual product.